

Particle Shadow Velocimetry System

User Manual

Table of Contents

PSV Background	.2
System Components	. 3
Network Setup	. 4
Experimental Setup	.7
Balancing the Signals	. 10
Data Acquisition	. 11
Post Processing	.13
Export Disclaimer	. 14

For questions or comments, please contact ISSI

Innovative Scientific Solutions, Incorporated 7610 McEwen Road Dayton, OH 45459 Ph.: (937) 630-3012 Fax: (937) 630-3015 **Tech Support:** <u>support@innssi.com</u> **Sales:** <u>issi-sales@innssi.com</u> **Website:** <u>www.innssi.com/psv</u>



Revision Date: 5/30/2019

PSV Background

Cost and safety issues associated with laser-based particle image velocimetry systems effectively eliminate their use in undergraduate student training. ISSI has developed a cost-effective alternative, Particle Shadow Velocimetry (PSV), illustrated below:



PSV uses light extinction (particle shadows) rather than light scattering in a laser sheet, resulting in high-quality particle imaging at much lower illumination levels. These safe illumination levels can be generated by a pulsed-LED system at a fraction of the cost of a pulsed laser. Multi-color LED illumination permits the use of a low-speed camera, further reducing system cost. Different color LEDs are sequentially pulsed, at high speed, and the sequential particle images are captured in a single frame on the camera's three color planes. Narrow depth-of-field optics are used to image a two-dimensional plane within the flow volume, similar to what is achieved with a laser sheet. Measurement close to surfaces is possible because the system is not degraded by surface reflections which severely limit laser-based systems. The PSV approach is fully compatible with the seeding approaches and specifications used with conventional PIV systems.



System Components

The PSV system is composed of a pulsed RGB LED, a CCD color camera, autofocus bellows, Canon® EF lens, Canon® EF lens controller and pulse delay generator.



Auto Focus Bellows

EF Zoom Lens

Network Setup

Begin by connecting each device to the Ethernet switch and connecting the power to each device. The pulse delay generator, camera and lens controller are all Ethernet controlled. Each should be directly connected via Ethernet to the switch. A fourth Ethernet cable should be connected to the computer so that all devices on the network can be accessed. Refer to the instruction manual for each device to set up communication and software. Next, connect BNC cables to the devices which will be triggered by the pulse delay generator. Connect channel 1 on the pulse delay generator to the camera trigger input. Connect channels 3 and 4 to the red and green BNC connections respectively on the LED module.



Network Configuration

The network on the PC NIC (network interface card) needs to be properly configured for communication over the network. To do this, navigate to the *'Network Connections'* page on the control panel where the local networks of the computer is displayed. Right-click on the network where the network switch is connected and select *'Properties.'*

From the 'Local Area Connection X Properties' window, click on 'Internet Protocol Version 4 (TCP/IPv4)' and then press the 'Properties' button, now useable. This will open the 'Internet Protocol Version 4 (TCP/IPv4) Properties' window. This is where the IP addresses are entered so that connected devices can be reached over the local network. The settings needed for communication are **IP Address** and **Subnet mask**.

Innovative Scientific Solutions, Inc.

The IP address of the NIC should use the following conventions:

IP Format:

Network.Network.Subnet.Host

- LC-2 IP: 192.168.2.252. Can also be set by user.
- PSG-3 IP: 192.168.2.30. Can also be set by user.
- Camera IP: Dynamic. Can be set by user.
- NIC IP: 192.168.2.XXX

TCP/IP Address	2 ×
IP address:	192 . 168 . 2 . 252
Subnet mask:	255.255.255.0
	Add Cancel

Local Area Connection 2 Properties	Internet Protocol Version 4 (TCP/IPv4) Properties
Networking	General
Connect using: Previous PCIe GBE Family Controller #2	You can get IP settings assigned automatically if your network supports this capability. Otherwise, you need to ask your network administrator for the appropriate IP settings.
Configure This connection uses the following items:	 Obtain an IP address automatically Use the following IP address:
Cleft of Microsoft Vetworks	IP address: 192 . 168 . 0 . 25
☑ ➡ QoS Packet Scheduler ☑ ➡ File and Printer Sharing for Microsoft Networks	Subnet mask: 255 . 255 . 255 . 0
	Default gateway:
A Internet Protocol Version 4 (TCP/IPv4)	Obtain DNS server address automatically O Use the following DNS server addresses:
Install Uninstall Properties	Preferred DNS server:
Description Allows your computer to access resources on a Microsoft network.	Alternate DNS server:
	Validate settings upon exit
OK Cancel	OK Cancel

The NIC IP address should have the same network and subnet addresses but a unique host, the last line of the IP address. The host can be any value 1-254 but different from the host address of the device. The subnet mask should be set to match that of the devices which, by default is 255.255.255.0. The **Default Gateway** and all other fields can be left blank.

Additional subnetworks can be configured for the same NIC. In the 'Internet Protocol Version 4 (TCP/IPv4)' window, the **Advanced** button will open the 'Advanced TCP/IP Settings' window. IP addresses of any subnetwork can be added for that NIC here. If multiple devices are being used on the same network and they have differing subnetworks, this option provides communication for all devices without the need to change IP addresses for each device.

If the lens controller is not connected properly a connection error (below) will be displayed. First, reset the device via the reset switch which will restores the default IP address.



If connection problems persist, check the connection of the Ethernet cable between the computer NIC and the each device on the network. To check that there is a physical connection, ping the IP address of each device from the Command Prompt. To do this, open the command prompt window and enter: **Ping XXX.XXX.XXX.-t**

C:\Windows\system32\cmd.exe - ping 192.168.2.109 -t	0 X	
Microsoft Windows [Version 6.1.7601] Copyright (c) 2009 Microsoft Corporation. All rights reserved.		*
C:\Users\Steve>ping 192.168.2.109 -t		-
Pinging 192.168.2.109 with 32 bytes of data: Reply from 192.168.2.109: bytes=32 time(ins TIL=128 Reply from 192.168.2.109: bytes=32 time(ins TIL=128		
		Ŧ

Experimental Setup

The camera and optics are placed opposite the pulsed LED with the flow field of interest in between. In PSV the setup permits the DOF, the field-of-view (FOV), and the working distance to be adjusted by introducing spacers or bellows between the camera body and lens. In a simple setup that uses bellows, the required extension of the bellows attachment is the product of the magnification and focal length.



The autofocus bellows allow for the Canon[®] EF lens controller to be used to control the focus and aperture of the attached Canon[®] EF lens. Adjustment of the bellows allows the field of view to be adjusted.



The pulse delay generator is used to control the pulse width, delay time and frequency of the camera and pulsed LED. A network switch is also included to communicate with each of the hardware components over the local network.



Camera Trigger Connection

Pulsed LED Trigger Connections

The pulsed LED is aligned 180-degrees from the camera lens. The LED is placed at a distance from the camera such that the LED is out of focus to the camera.



The LED must be out of focus relative to the camera so that the image background is near uniform. If the LED array is in focus relative to the camera, the background will be nonuniform and not provide the proper backdrop for particle shadow velocimetry. The LED and camera should be aligned to fill as much of the field of view as possible with an even background illumination.



LED Array Misaligned and in Focus

LED Array out of Focus

Above is an example of the correct focus and LED-camera distance so that the LED is completely out of focus.

Next, the flow field area should be focused. A printed target is used to focus the camera on the desired flow field location. Adjust the bellows on the lens so that the appropriate field of view comes into focus. Fine adjustments to focus can then be made by using the lens controller software. Once the field of view is in focus, the focus setting can be saved by the lens controller.



Target in Focus

Balancing the Signals

The signals from the red and green channels can be viewed in ProAcquire. Use the performance of the pulse width for each color output on the LED until the maximum signal is ~90% of dynamic range on each image.



Red Signal

Green Signal

Quick adjustments to the pulse width can be made by using the pulse delay generator software to adjust. The aperture can also be adjusted on the lens controller if needed. It is recommended the aperture be set fully open or near fully open to defocus the LED and minimize the depth of field.



LED Settings Lens

Controller Interface

For acquisition, the pulse delay generator is used as the master controller for the experiment. The camera should be set to external trigger, exposure control mode so that the timing settings are set by the pulse delay generator.



Channels 3 and 4 of the pulse delay generator will separately control the pulse width of the red and green channels and the delay time between them. The delay time between red and green will be dictated by the flow speed of the flow field studied.

Data Acquisition

Calibration Images

Prior to acquisition of data images, calibration images needed for post-processing should be acquired. Set the delay between red and green outputs on the RGB LED to 0.00. (On the pulse delay generator, set the delay time for the second pulse signal to the RGB LED to 0.00.) While running the seeded flow at a typical test condition, acquire ~50 images which will serve as calibration images. These will be post-processed using PIV software to create a color correction map, which will correct for pixel shift due to lateral chromatic aberration.



Zero Delay for Calibration Images



Calibration Image with Vectors Representing the Processed Correction Map

Flow Images

After calibration images are stored, images of the flow field can be acquired. The timing of the pulse delay generator will control the LED and camera acquisition of the flow field. The delay set between the red and green pulses of the LED should be such that an 8-10 pixel shift is observed between the red and green channels. This is dependent on the flow speed in the area of interest. Particle size should be 3-5 pixels



Red Signal

Green Signal

If the timing is set up correctly, the particle shift can be easily observed by taking a ratio of the red and green images.



Red-Green Shift

Post Processing

Commercial off-the-shelf PIV code can be used to process PSV data, after inverting the PSV images. First, the calibration images should be processed to get a mean vector field, which will likely radiate out from near the center of the image. This represents the chromatic aberration caused by the lens. After obtaining this field, the remaining images can be processed and the calibration vectors can be subtracted from the vectors in the data images to obtain the corrected vector field.

ISSI is developing a PSV Processing program for post-processing of PSV data, which will automatically perform this calibration. This program will be available in 2019 for the commercial PSV system.

Export Disclaimer

Any and all underlying information and technology contained in this document may be subject to U.S. export controls, including the Export Administration Act (50 U.S.C. Appx. §§ 2401 et seq.) and the Export Administration Regulations ("EAR", 50 C.F.R. Parts 730-774), and may be subject to export or import regulations in other countries. You are responsible for complying with all trade regulations and laws both foreign and domestic. Except as authorized by law or distributor agreement with ISSI, you agree and warrant not to export or re-export the information to any country, or to any person, entity, or end-user subject to U.S. export controls, including without limitation persons or entities listed on the U.S. Department of Commerce Bureau of Export Administration's Denied Parties List and the U.S. Department of Treasury's Specially Designated Nationals. You further represent and warrant that no U.S. federal agency has suspended, revoked, or denied your export privileges.